

Mycoplasma Removal Solution

Mycoplasma Removal Solution

Product Description:

CellBiologics Mycoplasma Removal Solution can be used to efficiently eliminate Mycoplasma contamination in the culture of mammalian cells. It can be used in combination with Antibiotic-Antimycotic Solution (Catalog No. PB-6920). Working concentration is 1:1000 dilution (Add 50µl Mycoplasma Removal Solution to 50 ml medium). Medium containing Mycoplasma Removal Solution and Antibiotic-Antimycotic Solution should be replaced every 2 days for two weeks.



Mycoplasma contaminations

Mycoplasma Removal Solution from *Cell Biologics* is distributed for in vitro research purposes only. Our products are not authorized for human use, for in vitro diagnostic procedures, or for therapeutic procedures.

Storage:

Store solution at 20°C.

How to use:

Add 50 µl Mycoplasma Removal Solution to 50 ml cell culture medium and replace the solution every second day for two weeks.

Product Information

Catalogue number: PB-6919 – 1 ml (sufficient for 5000 ml medium)
Species specificity: For Mycoplasma removal of mammalian cell cultures

ORDER UNDER sales@pelobiotech.com or FAX + 49 (0)89 517 286 59 88